FLEXCELL[®] FX-6000[™] Tension System

Stretch cells in monolayer or 3D culture

- Computer controlled bioreactor that applies **cyclic** or **static** tensile strains to cells cultured in monolayer or 3D allowing users to simulate *in vivo* tissue strains and frequencies.
- State-of-the-art digital valve automatically regulates and maintains regulated vacuum **and** positive air pressure for a specified regimen to deform a flexible-bottomed Flexcell[®] culture plate producing up to 33% substrate elongation (*see Table 1*).
- FlexSoft® software allows for multiple frequency, amplitude, and waveform changes to be programmed in one regimen.
- Cylindrical or Arctangle[®] loading posts used to apply equibiaxial or uniaxial strain, respectively.
- Gradient strain (unconstrained distention) achieved by removing the loading stations.
- Up to four FlexLink[®] controllers can be driven with one FX-6000[™] System.
- Available waveforms: static, sinusoidal, heart stimulation (Fig. 3), triangular, square, and custom.



Figure 3. FlexSoft® waveform plot showing a heart waveform.

Table 1.	Maximum achievable strains in Flexcell®
culture p	lates with FX -6000 $^{\text{\tiny TM}}$ Tension System.

Culture Plate	Max Strain
BioFlex® (with 25 mm Loading Stations)	21.8%
BioFlex® (without Loading Stations)	33%
Tissue Train®	20.8%
UniFlex®	12.2%
HT BioFlex®	8.0%

FX-6000[™] **Tension System includes:**

Equibiaxial Configuration

- Laptop with 15" monitor
- FlexSoft[®] FX-6000[™] software
- FX-6000[™] Tension FlexLink[®]
- BioFlex[®] Baseplate Kit
- Tubing and Adaptor Kit

Uniaxial Configuration*

- Laptop with 15" monitor
- FlexSoft® FX-6000[™] software
- FX-6000[™] Tension FlexLink[®]
- BioFlex® Baseplate Kit
- Tubing and Adaptor Kit
- TT4000 Set of Arctangle® Loading Stations
- *All components listed not included in price below

ORDERING INFORMATION (Please contact distributor for pricing)				
Catalog Number	Product/Item	Price		
FX-6000T	Flexcell® FX-6000™ Tension System			
FX-6000TFL	Additional FX-6000™ Tension FlexLink®			