4400-DOC001

System User Manual

Cryostation

CryoAdvance-50 | CryoAdvance-100

Version: 1.1 September 2022

www.montanainstruments.com support@montanainstruments.com





This page intentionally left blank.

Specifications and product information listed in this document are accurate at the time of publishing for a standard system. Options, custom designs, and other modifications may cause slight differences. Future design changes to the system, including software updates, may change information.

© Montana Instruments® Corporation, All Rights Reserved

No part of this publication may be reproduced, distributed, or transmitted in any form or by any means, including photocopying, recording, or other electronic or mechanical methods, without the prior written permission of Montana Instruments.

Montana Instruments, CryoAdvance, Cryostation, CryoOptic, and ATSM are trademarks of Montana Instruments Corporation. Other brand names used are the property of their respective owners.

Table of Contents

Section 1	- Preface	7
1.1	Conventions Used in this Manual	7
1.1.1	Abbreviations	7
1.1.2	International System of Units (SI) Symbols	8
1.1.3	System of Imperial Units Symbols	8
1.1.4	Explanation of Safety Warnings	8
1.1.5	Graphical Symbols	9
1.2	General Hazard Information	9
1.3	echnical Support Information	10
1.3.1	Warranty & Repairs	10
1.3.2	Accessories & Replacement Parts	11
1.3.3	Contact Details	11
Section 2	- System Overview	12
2.1	Cryostation	12
2.1.1	Intended Use	12
2.1.2	Components	12
2.1.3	Technical Specifications	13
2.1.4	Safety Information	15
2.2	system Control Unit	16
2.2.1	Intended Use	16
2.2.2	Components	16
2.2.3	Technical Specifications	19
2.2.4	Safety Information	19
2.3	/acuum Control Unit	20
2.3.1	Intended Use	20
2.3.2	Components	20
2.3.3	Technical Specifications	22
2.3.4	Safety Information	22
Section 3	- Sample Chamber Configuration	24
3.1	Accessing the Sample Space	24
3.2	Sample Mounting	25
3.3	Mounting a Thermometer	26

3.	4	Windows	28
	3.4.1	Window Covers	28
	3.4.2	Window Replacement	28
3.	5	Sample Chamber Wiring	31
	3.5.1	Types of Wiring	31
	3.5.2	Thermal Lagging Techniques	31
	3.5.3	User Wiring Interfaces	31
	3.5.4	Wiring Schematics	32
3.	6	System Care	36
	3.6.1	When Working in the Sample Chamber	36
	3.6.2	When Not Using The Cryostation	36
	3.6.3	Standard Checks Before Every Cooldown	36
	3.6.4	Periodic Checks Every 6-12 Months	37
3.	7	Long-Term Maintenance	37
	3.7.1	20,000 Hour Maintenance	37
Sect	tion 4	- System Usage & Operation	38
4.	1	System Control Options	38
	4.1.1	User Interface	38
	4.1.2	Network Interface	40
4.	2	Primary Operations	41
	4.2.1	Turning on the System	41
	4.2.2	Controlling System Temperature	42
	4.2.3	Venting and Pulling Vacuum	43
4.	3	Moving the System	44
Sect	tion 5	- Diagnostics & Troubleshooting	45
5.	1	Expected System Performance	45
5.	2	System Diagnostics	45
	5.2.1	Performance Issues	45
	5.2.2	Power and Communication Issues	46
	5.2.3	Compressor Issues & Fault Messages	47
5.	3	System Checks	49
	5.3.1	Temperature Optimization	50
	5.3.2	Vibration Mitigation	52

System User Manual

	5.3.3	Vacuum Check	52
	5.3.4	Helium Check	53
Sect	tion 6	- Appendices	. 55
6.		Related Documentation	

Section 1 - Preface

MARNING

Read all instructions before using this product

All users must read and understand this manual and all other safety instructions before using the equipment. Retain these instructions for future reference.

This manual is intended for users of Montana Instruments products and systems described herein. Users include anyone who may physically interact with the system or peripheral equipment, including installation, setup, or configuration. This includes anyone who may operate system components via control panels, the supplied user interface, or remote interfaces.

This manual may be used by facilities personnel for determining infrastructure requirements in the room or building where the equipment will be installed.

This manual is meant to be referenced by authorized service personnel for important safety and hazard information and other product restrictions.

1.1 Conventions Used in this Manual

The following style conventions are used in this document:

- Vertical bar (|)
 - o Indicates alternative selections. The bar may be used in place of "and" or "or".
- Alphanumeric List (1., 2., 3...| a., b., c...)
 - Indicates instructions or actions which should be completed in a specific ordered sequence.
- Bulleted List (• | | -)
 - o Indicates instructions, commands, or additional information about an action.
 - o May alternatively be used for unordered lists of materials or additional reference notes.
- Courier Font
 - o Indicates a label or indicator on a physical product or part.
 - o Indicates a system output, such as a display reading.
 - o May also be used for URLs, file paths, file names, scripting language, prompts, or syntax.

1.1.1 Abbreviations

The following abbreviations may be used:

- ACM: Ancillary Control Module
- CAN: Controller Area Network

- DMM: Digital Multimeter
- HDMI: High-Definition Multimedia Interface
- MI: Montana InstrumentsPCB: Printed Circuit Board
- TCM: Temperature Control Module
- UI: User Interface
- UPS: Uninterruptible Power Supply
- USB: Universal Serial Bus
- VNC: Virtual Network Computing

1.1.2 International System of Units (SI) Symbols

- C: Celsius
- cm: Centimeter
- K: Kelvin
- kg: Kilogram
- m: Meter
- mK: Millikelvin
- MPa: Megapascal
- mTorr: Millitorr
- mW: Milliwatt
- s: Second

1.1.3 System of Imperial Units Symbols

- ft, ': Foot
- in, ": Inch

1.1.4 Explanation of Safety Warnings

Safety and hazard information includes terms, symbols, warnings, and instructions used in this manual or on the equipment to alert users of precautions in the care, use, and handling of the system. The following hazard levels and information are considered:

A DANGER

Serious personal injury

Imminent hazards which, if not avoided, will result in serious injury or death.

MARNING

Serious personal injury

Potential hazards which, if not avoided, could result in serious injury or death.

A CAUTION

Possible personal injury

Potential hazards which, if not avoided, could result in minor or moderate injury.

NOTICE

Command or Product Safety Notice

Potential hazards which, if not avoided, could result in product damage.

» NOTE

Points of particular interest for more efficient or convenient equipment operation; additional information or explanation.

1.1.5 Graphical Symbols

The following symbols may be used in diagrams, supporting text, and on physical parts:

	Hazard Alert: General Warning	4	Hazard Alert: High Voltage
**	Hazard Alert: Laser Radiation	нэті	HDMI port
器	CAN bus module		USB port

1.2 General Hazard Information

The following descriptions are of general hazards and unsafe practices that may result in product damage, severe injury, or death.

- The products, parts, and components in this manual are to be serviced by authorized Montana Instruments service representatives only. Failure to do so will void the warranty and may damage the product and/or create a safety hazard.
- Only use components provided for the intended purpose described herein.
- If the equipment or any component is used or modified in a manner not specified by the manufacturer, the protection provided by the equipment may be impaired.

The following hazards may be typical for this product:

WARNING

Risk of injury when lifting or moving system components

System components, including standalone equipment and installed assemblies, may be heavy.

- Use caution when lifting or moving equipment or assemblies. Ensure proper lifting principles are used to avoid injury.
- Equipment or assemblies >20 kg should always be lifted by two or more people or with a suitable lifting device.

WARNING

High voltage: danger of electric shock

Electric shocks and burns from capacitor discharge or power circuits could lead to serious injury or death.

- Before turning on any power supply, the ground prong of the power cord plug must be properly
 connected to the ground connector of the wall outlet. The wall outlet must have a third prong or
 must be properly connected to an adapter that complies with these safety requirements.
- Only use replacement power cords or power plugs with the same polarity and power rating as that of the original ones. Do NOT use inadequately rated cables.

If the equipment or the wall outlet is damaged, the protective grounding could be disconnected.

- Do NOT use damaged equipment until its safety has been verified by authorized personnel.
- Do NOT disconnect or tamper with the operation of the protective earth terminal inside or outside the apparatus.

NOTICE

Only clean exterior surfaces with acceptable fluids

- Only use deionized water, glass cleaner, or isopropyl alcohol to clean the exterior surfaces of any
 enclosure. Do NOT use any volatile chemicals other than isopropyl alcohol.
- Apply fluid to a clean, lint-free cloth and wipe the surface with a cloth. Do NOT apply fluid directly to any surfaces or enclosures.

1.3 Technical Support Information

Any technical questions or issues with the system that cannot be resolved with the information in this manual should be referred to an authorized Montana Instruments service representative.

1.3.1 Warranty & Repairs

If the system or parts need to be returned to the Montana Instruments factory or an authorized service center for repair or service, contact an authorized service representative for a return merchandise authorization (RMA) number and instructions on returning the unit.

For a copy of the Limited Warranty Agreement, visit:

https://www.montanainstruments.com/support/warranty-information

1.3.2 Accessories & Replacement Parts

Only use cables, hoses, accessories, and parts provided or approved by the manufacturer. Follow all instructions for proper installation or replacement.

- To order spare or replacement parts, please contact your local service representative.
- To order new accessories or options, or for more information on other Montana Instruments products and technologies, please contact your local sales representative.

1.3.3 Contact Details

For a complete list of sales and service centers visit: www.montanainstruments.com/Contact

North American Authorized Service

- M-F 8:30am-5pm MST | Call: +1.406.551.2796
- Email: support@montanainstruments.com

North American Sales


- M-F 8:30am-5pm MST | Call: +1.406.551.2796
- Email: sales@montanainstruments.com

International Sales & Authorized Service

 Visit <u>www.montanainstruments.com/Contact</u> for contact information for your local representative.

Section 2 - System Overview

2.1 Cryostation

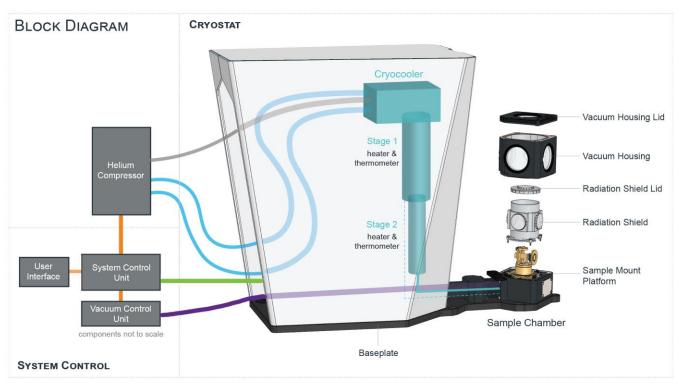

2.1.1 Intended Use

The system is a closed-cycle optical Cryostation used to control the temperature and vacuum parameters of an integrated sample environment.

The Cryostation is designed to connect to the system control unit, vacuum control unit, and helium compressor provided by Montana Instruments. The back panel of the Cryostation includes interface locations for these cable and hose connections.

2.1.2 Components

Cryostation

The Cryostation consists of the cooling tower assembly and the sample chamber, connected by a baseplate and a semi-rigid bellows assembly. The entire Cryostation assembly is rigidly mounted at either 45° or 90° to the hole pattern on an optical table via screw locations in the baseplate (optical table not included). The Cryostation can be mounted on either imperial or metric optical tables.

Cooling Tower Assembly

The main cooling is provided by a two-stage Gifford-McMahon (GM) cryocooler suspended inside the cooling tower assembly (the vertical aluminum cylinder inside the sheet-metal rear case). A vibration

damping support structure serves to isolate the cryocooler mechanical vibrations from the optical table and sample platform. The cryocooler is part of a closed-loop flow of helium which is pressurized by the separate helium compressor.

The cryocooler has two principal stages, each with thermometers to monitor temperature and heaters for warmup. Stage 1 is thermally coupled to the radiation shield inside the sample chamber. Stage 2 is thermally coupled to the sample mount platform.

Sample Chamber

The sample chamber consists of the lower vacuum housing with interfacing side panels, the sample mount platform and sample mount assembly, a surrounding radiation shield with inner "cold" windows, and the upper vacuum housing assembly with outer "warm" windows.

Thermal fluctuations are damped using both active and passive techniques. The sample platform mitigates the effects of gross thermal contraction during cooldown by using a thermal contraction-canceling design of the sample support structure.

Helium Compressor

Compressor unpacking and installation instructions are contained in the <u>Installation Procedure</u>. Refer to the Operating Instructions provided by the original manufacturer for further technical details, advanced operating instructions, and important safety information regarding the compressor. Pressure sensors allow for system monitoring and can aid in the detection of leaks.

Temperature Control Module (TCM)

A TCM peripheral card in the attached system control unit is used to read the thermometers and control the heaters connected to the cryocooler stages, platform, and user channels and feed that data to the appropriate readouts in the UI. This communication interface provides closed-loop control of the platform temperature.

The DSUB37 port (HEATER/THERMOMETER CONTROL) on the TCM card is used to interface with the Cryostation.



2.1.3 Technical Specifications

Environmental Specifications

Temperature of Environment	5 – 40 °C	
Humidity	5 – 80% non-condensing	
Altitude	<2000 m	

Power Specifications

	Model	Compressor: B-01
--	-------	------------------

Mains Power Con	nector on Unit	IEC 60320 C20
Line Voltage		208 – 240 VAC
Frequency		50 – 60 Hz
Maximum Current	t Draw	15 A
Maximum Power	Consumption	3.12 kW
Wall Outlet /	N. America & non-EU	NEMA 6-20R single-phase (see note below)
Receptacle	CEE Europe (non-UK)	CEE 7/3 or CEE 7/5 w/ common ground terminal
	UK	IEC 60309-6H (16 A)
	Israel	SI 32 (IS1-16P)

» NOTE

The NEMA 6-20R is a special outlet in the United States. An electrician must set up this circuit for single-phase 208-240 VAC with a 20 A breaker. **The circuit cannot be three-phase**.

Physical Dimensions

Component	L x W x H	Mass
Cryostation		
CryoAdvance-50	56.3 cm x 23.2 cm x 49 cm	29 kg
CryoAdvance-100	64 cm x 23.2 cm x 49 cm	29 kg
Compressor	61 cm x 38 cm x 64 cm	93 kg

2.1.4 Safety Information

The following hazards may be typical for this product:

MARNING

Risk of serious injury due to hazards associated with cryocooling

All personnel working with the system must be aware of the potential hazards associated with cryocooling.

• Personnel working with the system should be trained in emergency measures that may be required in the event of an accident.

Risk of suffocation due to potential asphyxiates.

Nitrogen (N_2) and Helium (He) are potential asphyxiates if released into an enclosed area with poor ventilation. A decrease in air oxygen content can be caused by leaks.

• Ensure that proper tubing is used, and good connections are made at each connection point to prevent the release of these gases.

Risk of explosion due to high pressure if the system is not allowed to vent properly.

• Never bolt or otherwise fasten the lid of the sample chamber closed. The lid acts as a safety pressure release in the event of high-pressure accumulation in the Cryostation.

Risk of cold contact burns.

Parts of this system are very cold and may cause severe burns to the skin.

Allow components to warm up to room temperature before touching. If contact occurs, consult a
physician immediately.

NOTICE

Take care when moving the Cryostation

- Do NOT tilt the Cryostation more than 45 degrees. Inverting the Cryostation will cause damage.
- The Cryostation and sample chamber are a single unit. The attached sample chamber must be supported at all times. Do NOT lift the Cryostation by the sample chamber.
- Do NOT lift the Cryostation by the cryocooler tube or the top of the main body enclosure.
- The Cryostation ships with red locking plugs and shipping support to prevent damage to sensitive components. Do NOT remove these until after the unit has been attached to the table.

Take care when moving the compressor

- Do NOT tilt the compressor. Doing so may damage the unit.
- The compressor is on casters for moving. Ensure casters are locked before operating.

Risk of product damage due to improper use

- Never disconnect the vacuum hose while the temperature of any stage of the Cryostation is below 285 K.
- Never open the case or vent valves when the temperature of any stage is below 285 K.
- Only use dry, high-purity nitrogen gas with the Cryostation. Do NOT substitute other gases for system venting.
- Avoid using any material in the sample chamber that may outgas or otherwise contaminate the optical windows and Cryostation surfaces.
- When manually operating heaters, monitor the Stage 1 and Stage 2 temperatures to ensure these temperatures do NOT rise above 350 K. Temperatures above 350 K may damage critical components within the system.

2.2 System Control Unit

Models SC1160





System Control Unit with User Interface Touchscreen

Front of Enclosure

2.2.1 Intended Use

The system control unit is a device used for instrumentation control of Montana Instruments products. It provides both the electronics hardware and software interface for communicating with other devices and controlling various instrument parameters.


2.2.2 Components

The system control unit consists of an outer enclosure, single-board computer, peripheral cards, a touchscreen user interface display, and several communication ports. The enclosure is compatible with a standard 19-inch-wide equipment rack (4U) and can be mounted with other rack-mount devices. It can alternatively be placed independently on a table or shelf.

User Interface Touchscreen

A 10-inch touchscreen display provides the main user interface control for the system. The software can alternatively be monitored and controlled via a VNC interface or remote scripting.

Enclosure Communication Ports

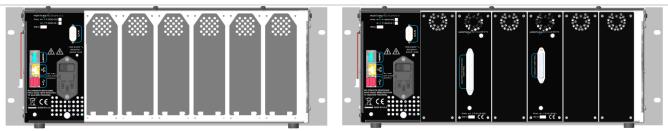

SC1160 (no peripheral cards) rear enclosure

SC1160 (two peripheral cards) rear enclosure

HDMI

The HDMI port on the rear face of the enclosure is used to interface with the touchscreen display.

USB

Two USB ports are available on the rear face of the enclosure. One of the USB ports is used to interface with the touchscreen display. The USB port is used to interface and communicate with the VC1130 turbopump option, if applicable.

Four USB ports are available on the front face of the enclosure. These can be used for connecting with external storage drives.

Ethernet

An ethernet port is available on the rear face of the enclosure for connecting with a local network or computer. A local network connection is required for controlling the instrument via the remote graphical display or scripting features, and a network internet connection is required for remote technical support.

DSUB9 (CAN)

The DSUB9 port on the rear face of the enclosure is used to interface with the helium compressor. This connection allows the system control unit to run the CAN bus in the compressor and display compressor readouts in the UI.

Peripheral Cards

The SC1160 contains slots for up to six, rear-facing peripheral control cards. There are several types of peripheral cards for controlling various equipment. The two standard cards are described below.

Ancillary Control Module (ACM)

An ACM peripheral card in the attached system control unit is used to control the valves and pump in the vacuum control unit. The DSUB25 port (VACUUM CONTROL) on the ACM is used to interface with the vacuum control unit.



Temperature Control Module (TCM)

A TCM peripheral card in the system control unit is the communication interface used to control the thermometry and heaters located on the cryocooler stages, platform, and user channels and display those readouts in the UI.

The DSUB37 port (HEATER/THERMOMETER CONTROL) on the TCM card connects to the DSUB37 port on the back panel of the Cryostation.



2.2.3 Technical Specifications

Environmental Specifications

Temperature of Environment	5 – 25 °C	
Humidity	5 – 80% non-condensing	
Elevation	<2000 m	

Power Specifications

	Model	SC1160	
Mains Power Conn	nector on Unit	IEC 60320 C14	
Line Voltage		100 – 240 VAC	
Frequency		50 – 60 Hz	
Maximum Current	Draw	6.65 A	
Maximum Power Consumption		665 W	
Wall Outlet /	/all Outlet / N. America & non-EU Standard NEMA 5-15		
Receptacle CEE Europe (non-UK)		CEE 7/3 or CEE 7/5 w/ common ground terminal	
	UK	BS1363 (UK) w/ common earth ground terminal	
	Israel	I-32-3 w/ common earth ground terminal	
	Cart Power Module	IEC 60320 C13	

Physical Dimensions

Component	LxWxH	Mass
SC1160	43 cm x 38 cm x 17 cm	13.6 kg

2.2.4 Safety Information

The following hazards may be typical for this product:

! WARNING

Risk of injury due to sharp edges

The interior of the enclosure contains sheet metal parts that may have sharp edges.

• When working inside the enclosure (authorized service personnel only), exercise caution to avoid injury.

MARNING

High voltage: danger of electric shock

Electric shocks and burns from capacitor discharge or power circuits could lead to serious injury or death.

- before accessing the enclosure or when otherwise servicing the unit (authorized service personnel only), completely power down the system and unplug the power cable.
- If power must be applied to diagnose issues or otherwise, a grounding strap must be applied to the arm interfacing internal components.

NOTICE

Peripheral cards must not exceed 600 W to avoid product damage

• The system control unit can supply a maximum power of 600 W across all installed peripheral cards. Ensure the cumulative power of all installed peripheral cards (maximum power rating of all cards added together) does not exceed 600 W.

Transportation and installation


- When not mounted in a rack unit, the enclosure should not be stacked on any other equipment, nor should other equipment be placed on it.
- Allow 8 cm minimum clearance from any ventilated face (sides, front) and 20 cm clearance in the rear for cables and hoses.
- Do NOT move the unit during operation. Disconnect all cables before moving. Lift the enclosure by using both handles on the front face.

2.3 Vacuum Control Unit

	Models	
VC1110		
VC1130		



Vacuum Control Unit

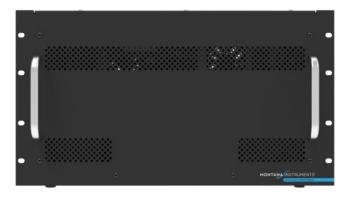

Front of Enclosure

2.3.1 Intended Use

The vacuum control unit is used for vacuum control of Montana Instruments Cryostations. It is designed to connect to the system control unit and Cryostation provided by Montana Instruments. It is recommended to locate the vacuum control unit near both connected apparatuses.

2.3.2 Components

The vacuum control unit consists of an outer enclosure, diaphragm (roughing) pump, ¼ inch vacuum coupling O-ring (VCO) connection, one case valve, one vent valve, and several communication ports. The enclosure (6U) is compatible with a standard 19-inch-wide rack and can be mounted with other rack-mount devices. It can alternatively be placed independently on a table or shelf.

The roughing pump is used to achieve a low vacuum. A high vacuum is achieved with cryopumping. During cooldown, the sample platform temperature lags the Stage 1 temperature in the Cryostation by ~100 K this allows for remaining molecules to be trapped by the charcoal adsorbers and cold surfaces conductively connected to Stage 1 rather than the sample. This minimizes sample contamination.

Enclosure Communication Ports



VC1110 rear enclosure

VC1130 (Turbo) rear enclosure

DSUB25 (VACUUM CONTROL)

The DSUB25 port on the rear face of the enclosure is used to interface with the ACM in the system control unit. This connection allows the system control unit to control the pumps and valves in the vacuum control unit and display the status of the roughing pump, vent valve, case valve, and system pressure readouts in the UI.

USB (VC1130 only)

The USB port on the rear face of the enclosure (VC1130 turbopump version only) is used to interface with the system control unit, allowing the turbopump to be controlled via the UI.

Ethernet (VC1130 only)

The ethernet port on the rear face of the enclosure (VC1130 turbopump version only) is used for communicating with the turbopump vacuum gauge.

2.3.3 Technical Specifications

Environmental Specifications

Temperature of Environment	5 – 25 °C
Humidity	5 – 80% non-condensing
Elevation	<2000 m

Power Specifications

	Model	VC1110, VC1130
Mains Power Conr	nector on Unit	IEC 60320 C14
Line Voltage		100 – 240 VAC
Frequency		50 – 60 Hz
Maximum Current	Draw	1.9 A
Maximum Power C	Consumption	190 W
Wall Outlet /	N. America & non-EU	Standard NEMA 5-15
Receptacle	CEE Europe (non-UK)	CEE 7/3 or CEE 7/5 w/ common ground terminal
	UK	BS1363 (UK) w/ common earth ground terminal
	Israel	I-32-3 w/ common earth ground terminal
	Cart Power Module	IEC 60320 C13

Physical Dimensions

Component	LxWxH	Mass
VC1110 VC1130	43 cm x 38 cm x 26.5 cm	24 kg

2.3.4 Safety Information

The following hazards may be typical for this product:

MARNING

Refer to associated product manuals for complete safety information

• The VC1130 model contains a Pfeiffer Vacuum® HiPace® 80 turbopump. Refer to the Operating Instructions provided by Pfeiffer for important safety information regarding this component.

MARNING

Risk of injury due to sharp edges

The interior of the enclosure contains sheet metal parts that may have sharp edges.

 When working inside the enclosure (authorized service personnel only), exercise caution to avoid injury.

WARNING

High voltage: danger of electric shock

Electric shocks and burns from capacitor discharge or power circuits could lead to serious injury or death.

- Completely power down the system and unplug the power cable before accessing the enclosure or when otherwise servicing the unit (authorized service personnel only),.
- If power must be applied to diagnose issues or otherwise, a grounding strap must be applied to the arm interfacing internal components.

NOTICE

Risk of product damage due to improper use

- Do NOT kink vacuum tubing or install vacuum tubing in an area where it may be pinched.
- Ensure all vacuum clamps and fittings are secured tightly before operation.
- Do NOT disturb vacuum tubing while the turbopump (if applicable) is operational.

Risk of turbopump damage due to improper use (VC1130 model only)

Venting the system at rotor speeds higher than 250 Hz may cause damage to the turbopump

- Do NOT vent the system when the turbopump is rotating faster than 250 Hz. Observe rotational speed on the UI to ensure safe rotating speeds before manual venting.
- If the system control unit unexpectedly shuts down before turning off the turbopump, turn off the main power to the vacuum control unit and wait at least 15 minutes before venting the system.

Transportation and installation

- When not in a rack unit, the enclosure should not be stacked on any other equipment, nor should other equipment be placed on it.
- Allow 8 cm minimum clearance from any ventilated face (sides, front) and 20 cm clearance in the rear for cables and hoses. For the VC1130, allow 30 cm rear clearance.
- Do NOT move the unit during operation. Disconnect all cables before before moving. Lift the
 enclosure by using both handles on the front face.

Section 3 - Sample Chamber Configuration

3.1 Accessing the Sample Space

Before starting the Cryostation, check the sample space. Depending on the configuration, there may be foam inserts to remove before cooling the system down for the first time. Check the Manual Addendum for instructions specific to your configuration.

NOTICE

Keep sample chamber and surfaces clean


- Always wear clean gloves when working in the sample chamber to avoid contaminating surfaces.
- Take care not to touch the optical windows on the lid, vacuum housing, or the internal radiation shield. Window covers are provided to keep the windows clean.
- Keep O-ring seals clean and free from debris. Do NOT place the housing with an o-ring or sealing surface positioned downwards unless there are protruding bosses to keep it from touching the surface.
- When working inside the sample chamber, use extreme caution not to drop screws into the chamber assembly.

Start with the system at room temperature and the chamber vented to atmosphere.

- 1. Carefully lift the top lid from the sample chamber and place it on the table with the protruding bosses down. The lid may sit on the bosses without compromising the clean surface inside the lid.
- 2. Using two hands, carefully lift the window assembly from the sample chamber. For CryoAdvance-50: place the window assembly on the table so that it rests on a flattened corner (not on a window), this will reduce contamination and keep the windows free from scratches. The CryoAdvance-100 model includes bosses around the windows of the vacuum housing so that it may be placed on the table with the protruding bosses down.





Step 1: Lifting off the lid with protruding bosses on the underside

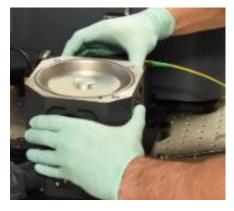


Step 2: Lifting off the window assembly

- 3. Lift the cap from the radiation shield to remove it.
- 4. Unfasten the 4x M3 button head screws on each side of the radiation shield; lift up and over any sample mount structures to remove.

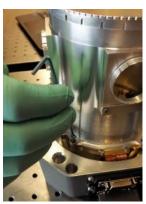

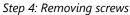

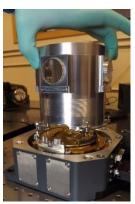

Step 4: Lifting off radiation shield assembly

NOTICE

Take care when reassembling the sample chamber

- Before replacing the vacuum housing & lid, carefully check O-rings and sealing surfaces for any loose fibers or debris (contamination could affect vacuum performance).
- Take care not to overtighten screws when replacing the radiation shield. Use of a torque screwdriver set to 5 in-lb is recommended.

3.2 Sample Mounting

Most systems ship with a user-specified sample mount, depending on the option(s) purchased.

- 1. Remove the previous sample from the face of the mount if needed. Clean off residual grease with a Kimwipe, lens tissue, or lint-free cloth (use a small amount of isopropanol or acetone if needed)
- 2. Apply a thin layer of new N-grease to the sample mount surface. Use a flat tool (not metal or cotton) to spread the grease out evenly, taking care not to scratch the surface.

NOTICE


Do not use N-grease for temperatures above 300 K

Apiezon® N-grease softens at 305 K. If the user plans to operate above 300 K, it is recommended to use an alternative sample mounting material to ensure a vertically mounted sample does not slide off the sample mount.


» NOTE

Depending on the sample, VGE, silver paint/paste, or copper SEM tape can be used in place of N-grease.


3. Using tweezers, set your sample on top of the grease layer and press down gently to ensure a good thermal connection.

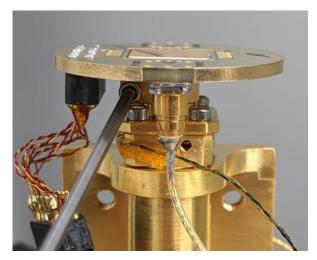


Attach sample mount to exchange boss

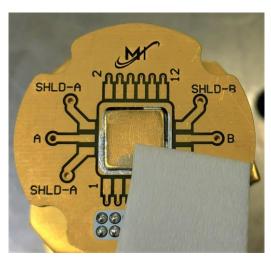

Step 1: Clean the surface of the mounting platform

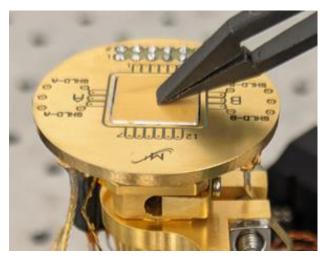

Step 2: Grease mounting surface

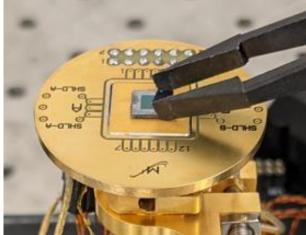

Step 3: Attaching new sample

» NOTE

To check if the sample is secure, hold the sample mount horizontally so the sample faces down and forcibly tap the mount with a finger.

3.3 Mounting a Thermometer

Many users choose to remove the sample thermometer and mount it elsewhere in the system or install a user thermometer. To do this:

NOTICE

Cernox® thermometer wires are extremely fragile.

Handle wires with care to avoid breaking.

1. Remove any VGE (the adhesive lagging the thermometer wire) with isopropyl alcohol or acetone. Clean off any residual VGE.

- 2. Clean off the area where the thermometer will be mounted with isopropyl alcohol or acetone.
- 3. Attach the thermometer, ensuring the bottom is in even, flat contact with the mounting surface.
 - a. **With Clamp:** Apply a thin layer of N-grease to the bottom of the thermometer. Clamp it down securely, ensuring no part of the thermometer "springs up" from the surface.
 - b. **Without Clamp:** Apply a small amount of VGE (thinned with acetone or isopropyl alcohol) to the bottom of the thermometer, wait 10 seconds, then place on the mounting surface. Use a small weight to apply pressure and hold the sensor in place while it cures (24 hours under ambient conditions or 30 minutes under 60 °C heat).



Step 3a: Thermometer clamped to the mounting surface

Wire lagged with VGE (clamp style)




Step 3a: Thermometer clamped to the mounting surface



Wire lagged with VGE (plug style)



Step 3c: Thermometer installed in a plug

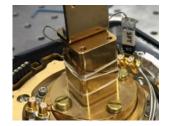

Wire tied in place with floss (plug style)

4. Tie the thermometer wire in place using unwaxed floss, nylon string, or Kapton® tape (optional). Apply VGE to approximately the first 0.5 in. (1 cm) of wire and press to a metal surface to properly lag the wire. Do not encase the wire in VGE.

» NOTE

- Be sure the thermometer and wire are completely dry before pulling vacuum or cooling down the system. The wire should not touch the radiation shield or sample mount.
- The sample temperature may read up to 0.2 K higher than the platform.

3.4 Windows

3.4.1 Window Covers

The standard sample chamber has five optical access locations. The system ships with window covers in place. The window covers are removed by turning counter-clockwise until the notches are aligned with the openings in the window retaining ring. To replace the covers, re-align the notches and turn clockwise until the cover locks in place.



» NOTE


Use care when removing the window covers if the system is running, as this may loosen the retaining ring holding the window in place.

3.4.2 Window Replacement


Vacuum Windows

For replacing outer (warm) vacuum windows, a special window tool is provided in the accessory kit:


- 1. Align the tabs on the window tool with the corresponding notches in the Delrin retaining ring holding the window in place.
- 2. Use the window tool to loosen the retaining ring by turning counter-clockwise.
- 3. Remove the retaining ring to access the window.
- 4. before reinstalling the window, check the O-ring to ensure it is clean and free of debris and has a very thin layer of L-grease (just enough so the surface is shiny).
- 5. Carefully center the window over the O-ring during reinstallation. Ensure that the O-ring is fully covered by the window. Avoid repositioning the window, as this may spread L-grease onto the clear aperture

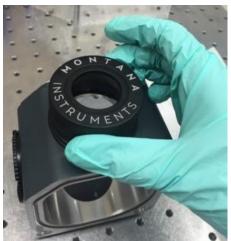

Steps 1-2: Using window tool to unscrew the retaining ring

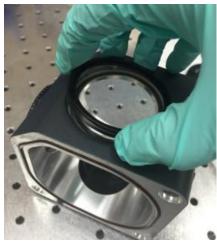

Step 3: Removing the retaining ring

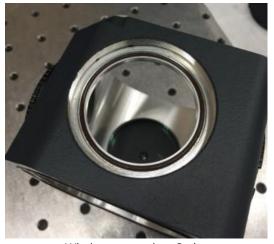

Window centered on O-ring

NOTICE


Take care when handling and removing windows

- Handle the radiation shield and windows carefully to prevent scratches or fingerprints.
- Before using the window tool, remove the housing from the system.
- Do not use too much force on the window tool as this could cause it to slip and damage the window.

Radiation Shield Windows

The inner (cold) windows on the radiation shield are held in place with a tension ring.

- 1. Depending on the model, the radiation window holders are removed by unscrewing the threaded tension ring or removing the four M2 screws surrounding the window.
- 2. Remove the window from the tension ring by pressing it out of the spring fingers. Some force may be required.
- 3. before reinstalling the window, add a very thin layer of N-grease to the edge of the spring fingers. This layer should be thin enough so none of the grease spreads onto the optic when reassembled.
- 4. Reinstall the window by press-fitting the tension ring onto the window. Ensure the spring fingers hold the window tightly and evenly.
- 5. For threaded tension rings, add a small amount of N-grease to the threads before screwing back in place.

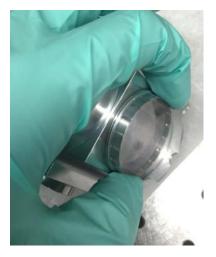

Step 1: Unscrewing threaded tension ring



Pressing window out of tension ring



Pressing tension ring onto the window

» NOTE

- The threaded holders can be adjusted to allow the sample to be positioned closer to external optics. Always ensure that no portion of the radiation shield touches the outer housing and that the radiation and vacuum windows do not touch.
- If a radiation window is removed, the added heat load can be as much as 50 100 mW per window. Depending upon the emissivity and thermal conductivity of the sample, local heating of the sample from this excessive radiation can cause the sample to be significantly warmer than the cold platform it is mounted to.

3.5 Sample Chamber Wiring

3.5.1 Types of Wiring

Various applications require different types of wiring. Always select the optimal wire for the application with the appropriate diameter (between 32-40 gauge). Longer wires are ideal -- the length of wiring between stages should be at least 2 in (5 cm).

- **32 AWG Phosphor Bronze wire**: low thermal conductivity and good for moderate power transmission
 - o Recommended use: Building additional wiring harnesses
 - o Raises temperature ~5.0 mK per wire
- Copper wire: high thermal conductivity and good for high power transmission
 - Recommended use: Wiring harness for high power transmission when high base temperatures are acceptable
 - o Raises temperature ~300 mK per wire

» NOTE

See our Cryostation Wiring Guide for details, theory, and illustrations.

3.5.2 Thermal Lagging Techniques

To minimize the effects of heat loads through wiring on the base temperature, all incoming wiring (including coax cables) must be properly thermally lagged to Stage 1 of the cryocooler via the radiation shield.

CryoAdvance-100 model provides "pre-lagged" connections for some user inputs and wiring. These wires, which plug directly into the 4 K circuit board "wedge" inside the sample space, do not need additional thermal lagging.

Wiring must route under the thermal clamp locations when pre-lagged connections are not available:

- 1. Unscrew the two M2 screws on the top of the thermal clamp. Remove the top cover.
- 2. Insulate the wires with a small piece of Kapton® tape to help prevent electrical shorts.
- 3. Place the wires on the felt pad. Wires should not cross or touch one another under the clamp.
- 4. Replace the top cover of the clamp and replace the two screws. Do NOT overtighten.
- 5. Check to ensure wires do not touch the inside of the radiation shield.

» NOTE

Ensure there are a few inches of wire before and after the thermal clamp. A 10-inch (25 cm) wire should be thermally lagged such that a 5.0 in. (12.5 cm) length is outside of the thermal clamp and the remaining 5.0 in. (12.5 cm) section is inside of the 4K sample space.

3.5.3 User Wiring Interfaces

The Cryostation provides integrated connectors for user DC feedthroughs, user thermometers, and heaters.

NOTICE

Do NOT remove sample chamber circuit boards

- The standard 4 K circuit board "wedge" (contains USER, SAMPLE THERM, etc.) used on CryoAdvance-100 models should not be removed as it can lead to the disconnecting of the platform thermostat and heater. If this wedge needs to be removed for any reason, please contact an authorized service representative for detailed instructions.
- The black printed circuit board on the CryoAdvance -50 model should not be removed. The black Delrin screws holding it in place should remain loose. Overtightening can cause the board to flex and make it difficult to obtain a proper vacuum seal.
- The Operational cable and associated wiring must remain connected for the system to be operational. Do NOT modify these connections.

User DC Connections

User DC connections allow users to route wiring to header pins on the sample chamber circuit board. These available connections can be used to interface with external connections and devices. The accessory kit comes with pin connectors for interfacing user inputs to the external connectors.

User Temperature Channels

User thermometer and heater connections allow users to add additional thermometers and heaters at header pin locations on the sample chamber circuit board. The system control unit provides open temperature channels (User 1 & 2) for manually controlling additional thermometers and heaters via the UI.

» NOTE


The provided sample thermometer is symmetric, so it can be connected to Pin 1 in either orientation.

3.5.4 Wiring Schematics

The wiring schematics for each model are outlined on the following pages.

CryoAdvance-50

All wiring must be routed from the sample chamber circuit board under the thermal clamps into the cold space.

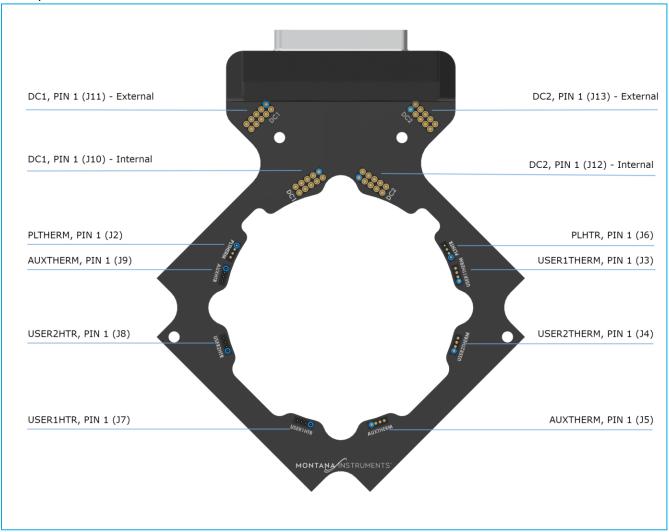


CryoAdvance-50 PCB diagram

User Interface	Internal Connections	
Temperature Channel	PCB Pins	
		PIN 1
		PIN 2
	USER1THERM	PIN 3
LICED 1		PIN 4
USER 1		PIN 1
	LICEDALITE	PIN 2
	USER1HTR	PIN 3
		PIN 4
		PIN 1
	USER2HTR	PIN 2
		PIN 3
LICED 3		PIN 4
USER 2	USER2HTR	PIN 1
		PIN 2
		PIN 3
		PIN 4
	AUXHTR	PIN 1
		PIN 2
		PIN 3
ALIV		PIN 4
AUX	AUXHTR	PIN 1
		PIN 2
		PIN 3
		PIN 4
User Temperature Channels		

30K Platform Internal	External	30K Platform Internal	External
DC1, PIN 1	DC1, PIN 1	DC2, PIN 1	DC2, PIN 1
DC1, PIN 2	DC1, PIN 2	DC2, PIN 2	DC2, PIN 2
DC1, PIN 3	DC1, PIN 3	DC2, PIN 3	DC2, PIN 3
DC1, PIN 4	DC1, PIN 4	DC2, PIN 4	DC2, PIN 4
DC1, PIN 5	DC1, PIN 5	DC2, PIN 5	DC2, PIN 5
DC1, PIN 6	DC1, PIN 6	DC2, PIN 6	DC2, PIN 6
DC1, PIN 7	DC1, PIN 7	DC2, PIN 7	DC2, PIN 7
DC1, PIN 8	DC1, PIN 8	DC2, PIN 8	DC2, PIN 8
DC1, PIN 9	DC1, PIN 9	DC2, PIN 9	DC2, PIN 9
DC1, PIN 10	DC1, PIN 10	DC2, PIN 10	DC2, PIN 10
User DC Connections			

CryoAdvance-100

Cryostation CA-100 user wiring to the 4 K circuit board shown below is "pre-lagged" and does not need to be routed under the thermal clamps. Other inputs will require proper thermal lagging.

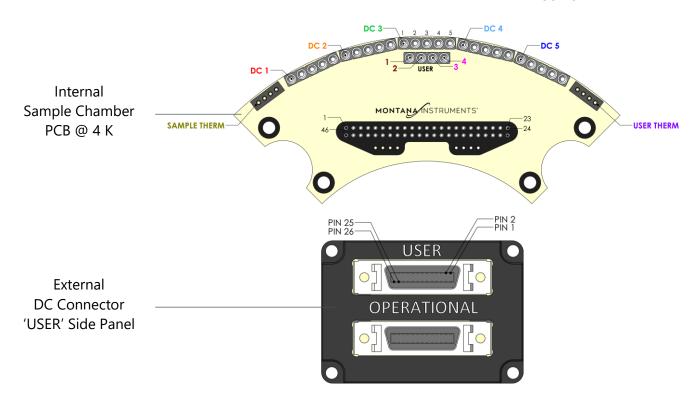

User Temperature Channels

User Interface		Internal Connections	
Temperature Channel		PCB	Pins
			PIN 1
	Thermo-	SAMPLE	PIN 2
USER 1	meter	THERM	PIN 3
USER I			PIN 4
	Haatau	LICED	PIN 3
	Heater	USER	PIN 4
	Haatau	LICED	PIN 1
USER 2*	Heater	USER	PIN 2
			PIN 1
-or- CRYO- OPTIC (if available)	Thermo-	USER	PIN 2
	meter	THERM	PIN 3
			PIN 4

^{*}With a Cryo-Optic installed, the USER 2 temperature channel and corresponding pins are not available to the user.

User DC Connections

PCB	Side Panel	РСВ	Side Panel
DC1, PIN 1	PIN 1	DC4, PIN 1	PIN 16
DC1, PIN 2	PIN 2	DC4, PIN 2	PIN 17
DC1, PIN 3	PIN 3	DC4, PIN 3	PIN 18
DC1, PIN 4	PIN 4	DC4, PIN 4	PIN 19
DC1, PIN 5	PIN 5	DC4, PIN 5	PIN 20
DC2, PIN 1	PIN 6	DC5, PIN 1	PIN 21
DC2, PIN 2	PIN 7	DC5, PIN 2	PIN 22
DC2, PIN 3	PIN 8	DC5, PIN 3	PIN 23
DC2, PIN 4	PIN 9	DC5, PIN 4	PIN 24
DC2, PIN 5	PIN 10	DC5, PIN 5	PIN 25
DC3, PIN 1	PIN 11	empty	PIN 26
DC3, PIN 2	PIN 12		
DC3, PIN 3	PIN 13		
DC3, PIN 4	PIN 14		
DC3, PIN 5	PIN 15		

3.6 System Care

Recommend system care procedures should be followed by any users of the system. For further information on any of these procedures, contact an authorized service representative for assistance.

3.6.1 When Working in the Sample Chamber

- Keep surfaces clean. Avoid touching any surfaces inside the sample space with your fingers as oils or other foreign contaminates can easily be transferred to the surfaces, the sample, or optics. **Always wear clean gloves.**
- Use proper grease and adhesives in the sample chamber. The accessory kit includes Apiezon® L-grease, N-grease, and the adhesive GE Varnish (VGE).
- Avoid using too much grease a thin layer (just enough so the surface is shiny) is best for metal-to-metal surfaces, samples, and O-rings. Excessive grease can increase outgassing and contamination of other surfaces in the sample chamber.
- Inspect, wipe, and grease O-rings. Make sure that the O-rings are clean with a thin layer of L-grease. The exposed surface should be wiped with a dry Kimwipe or lens tissue and re-greased every 10-15 uses.
- Check to ensure wires are preserved: make sure wires do not overlap under thermal clamps and that the clamps are not overtightened; make sure wires do not contact the radiation shield or the sample mount (aside from the wire bond pads).

3.6.2 When Not Using The Cryostation

- Keep the lid on the sample chamber to keep it free from dust.
- Supply a small amount of nitrogen to keep the system clean and dry.
 - Use the VENT command with "Vent Continuously" enabled to keep nitrogen flowing through the chamber at atmospheric conditions.
- Keep the sample chamber under vacuum. Use the PULL VACUUM command button to pull and keep the chamber under a medium vacuum state.

3.6.3 Standard Checks Before Every Cooldown

- Ensure the radiation shield is bolted down and secured. Make sure all radiation shield windows are in place.
- Ensure the sample is mounted properly.
- Ensure wiring from sample thermometers or other internal components does not touch the inside of the radiation shield.
- Check that the software starts up and reads all thermometers. The temperature channel values should be fluctuating a small amount, indicating the thermometer readings are active.

•	Check helium	pressure to e	ensure values	are within a	ppropriate \	values per	the table below:

Configuration	Equalization Pressure
F20-L, any hose length	1447 kPa (210 psig)
HC-4A2 with ½" dia. X ≤20 ft. flex hoses	1655 – 1670 kPa (240 – 245 psig)
HC-4A2 with ½" dia. X >20 ft. and ≤66 ft. flex hoses	1450 – 1480 kPa (210 – 215 psig)

3.6.4 Periodic Checks Every 6-12 Months

- Inspect VGE joints and repair if needed. VGE can flake off after several thermal cycles.
 - 1. Clean off old VGE with acetone or isopropyl alcohol then add a thin layer of VGE in the same location.
- Discard and replace VGE annually VGE has a limited shelf life of one year.

» NOTE

O-rings rarely need to be replaced unless they are nicked or damaged. If damaged, use Teflon-tipped tweezers or your fingers to remove the O-ring. Do NOT use metal-tipped tweezers or other sharp objects as this could damage the aluminum housing and create leaks.

3.7 Long-Term Maintenance

3.7.1 20,000 Hour Maintenance

The maintenance procedures below must be completed by a certified technician. This can be done onsite without removing the equipment or disturbing the optical setup. Contact an authorized service representative to schedule maintenance.

Cryocooler

The GM cryocooler normally needs maintenance every 20,000 hours of operation when driven by the variable speed compressor. After extended use, some of the bearings and bushings will wear down.

Other indicators of needing cryocooler maintenance include:

- The system has trouble cooling, or the achievable base temperature drastically degrades.
- The system makes a cogging or grinding noise in the cryocooler that persists continually.

» NOTE
Periodic chirping as the cryocooler runs is normal as the bearings wear in.

Compressor Filter and Adsorber


Refer to the system manual for the applicable SHI compressor model.

Section 4 - System Usage & Operation

4.1 System Control Options

4.1.1 User Interface

The touchscreen display provides the primary user interface control of the system.

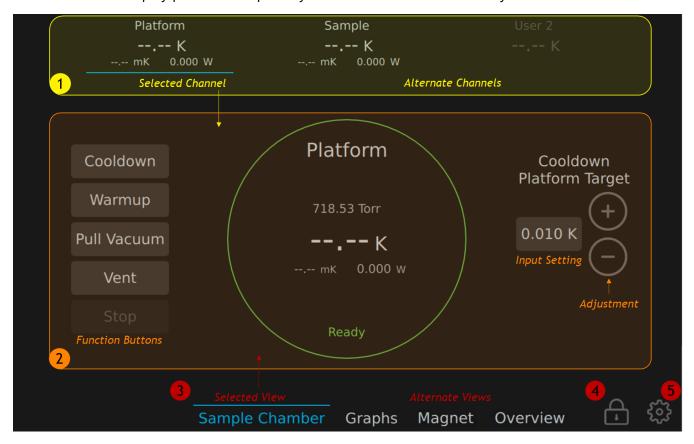

General Navigation

The example screen above shows the general layout, navigation, and controls for the UI. Not all controls and views will be available for all instruments.

- 1. **All Channels View:** Displays live status readouts of temperature, stability, and heater power for all attached channels of a given instrument. Press a channel to bring up its operation controls and settings. Channel names can be customized in the MENU.
- 2. **Selected Channel Operational View:** Displays the live status readouts and associated controls for the selected channel.
 - a. Function Buttons: Buttons with action statements tell the system to do something.
 - b. Adjustment Buttons: Circular buttons with icons provide operational control.

- c. **Input Settings:** Buttons with numerical values and units show the current input settings for a given command. To set new values, use the adjustment controls (if available) or press the display box to open an input popup dialog.
- 3. **Display Screens:** Choose which screen view to display. Alternate screens (when available) may show different readouts or have different control settings.
- 4. **Screen Lock:** Lock the touchscreen.
- 5. **Menu:** View or customize instrument and system settings. Other sub-menu features include:
 - a. **Remote Connections:** VNC and scripting parameters.
 - b. Data: Download system data onto a connected USB Flash Drive.
 - c. **Event Log:** Historical log of system-level events for reference or diagnostics.
 - d. **Maintenance:** Displays any recommended maintenance procedures.
 - e. **Tech Support:** Settings for assisting technical support personnel.

» NOTE

- Press and hold a command or input settings button to show on-screen help for that operation.
 Drag off the button before releasing to avoid executing the operation.
- The control buttons do not have on-screen help. Pressing and holding will execute the associated operation.

Application UI: Cryostation

Channel Types

- **Temperature Channel**: Displays live readouts of thermometers (temperature, temperature stability) and heaters (applied power) in the system.
 - Platform: The primary channel used to control the system. These commands will drive control operations across the Cryostation, compressor, and vacuum control unit.
 - User 1: The User 1 temperature channel displays the sample thermometer temperature readout.

» NOTE

- If the sample thermometer is not being used, it is recommended to turn it off to reduce noise on other temperature channels.
 - Navigate to MENU > INSTRUMENT SETTINGS > USER 1. Toggle the Temperature Channel Enabled switch to the left to disable.

Display Screens

- **Sample Chamber:** Displays temperature channel controls and readouts. This screen is used to run most primary system operations.
- Graphs: Displays real-time system data in graphical form.
- **Overview:** Displays system status readouts for all connected sub-systems. Press any value to pull up detailed information and additional control settings (if available) for that parameter.
 - o **Vacuum System:** status of pumps | valves | N₂ input | pressure gauges
 - o **Cooler:** status of the cryocooler/compressor | Stage 1 and Stage 2 temperature channels
 - o **Sample Chamber:** status of all sample chamber temperature channels

4.1.2 Network Interface

To utilize any of the provided network services, the system control unit must be connected to a local network using the Ethernet port on the rear face of the enclosure.

Each service uses a different network port number for communication. You may need to work with your IT department to configure these ports to be allowed through the firewall on your network.

To find your system's IP or MAC address:

- 1. On the touchscreen UI, navigate to MENU > REMOTE CONNECTIONS
- 2. Scroll down to find the system's unique IP and MAC addresses. Your IT department may need to configure the network to provide network access to the device.

Webserver | Onboard Documentation

♣ Standard HTTP port 80

The onboard webserver provides direct access to system-specific documentation, instrument scripting documentation and examples, web-based data dashboard links, select service documents, and other support materials.

To access the onboard documentation:

1. From an external computer connected to the same local network, use a web browser to navigate to the system's IP address (example: http://102.123.1.34).

Remote Graphical Display | VNC Viewer

- Non-standard HTTP port 2224

The user interface can alternatively be viewed or controlled via an external computer using Virtual Network Computing (VNC) technology. The external computer must have a VNC viewer program installed (several free options are available, including www.realvnc.com and www.tightvnc.com). When Remote Graphical Display is enabled, the VNC connection will mirror the UI control screen.

To enable Remote Graphical Display:

- 1. On the touchscreen UI, navigate to MENU > REMOTE CONNECTIONS
- 2. Toggle the Remote Graphical Display switch to ENABLED
- 3. Follow the onscreen directions for connecting to the IP address using a VNC viewer software

The VNC connection is password protected using a unique password for each system. The password is provided with your system documentation and in the onboard file QUICK START GUIDE - NETWORK INTERFACE CONNECTIONS. If lost, contact an authorized service representative to retrieve a copy.

Remote Scripting

- Non-standard HTTP port 47101

The instrument can alternatively be programmatically controlled via scripting commands using an external computer.

To use external scripting with the instrument:

1. On the touchscreen UI, navigate to MENU > REMOTE CONNECTIONS

- 2. Toggle the Instrument Scripting switch for the desired instrument to ENABLED
- 3. Follow the onscreen directions for opening the onboard scripting documentation, then use the available instructions for scripting with the instrument.

Remote scripting commands can be sent using an encrypted SSH tunnel. Refer to the onboard Instrument Scripting documentation on the webserver for further details.

Data Dashboards

♣ Non-standard HTTP port 8050

Data dashboards allow the user to view and analyze system data from a web browser.

To enable and view data dashboards:

- 1. On the touchscreen UI, navigate to MENU > REMOTE CONNECTIONS
- 2. Toggle the Data Dashboards switch to ENABLED
- 3. From the onboard documentation, click the SYSTEM DATA DASHBOARD link to access the system data.

MI Remote Support

- Secure SSH tunnel port 2223

The Montana Instruments Remote Support connection is used by authorized service representatives to remotely access your system to provide diagnostic and troubleshooting support. If the system is behind a NAT/Firewall, the firewall must allow the system to connect a Secure SSH tunnel on port 2223 to mirs.montanainstruments.com.

4.2 Primary Operations

4.2.1 Turning on the System

Before starting the power-up procedure, ensure that the power switches on the back of the system control unit and the front of the compressor are toggled to OFF (o).

- 1. If a dry nitrogen source is connected to the system, verify the supply pressure is set to ~15 psi.
- 2. Ensure the black circuit breaker switch on the front of the compressor unit is toggled to ON (|).

NOTICE

The green power switch should remain OFF (o) during operation as this is controlled by the remote ON/OFF control.

- 3. Turn on power to the vacuum control unit by toggling the power switch on the back of the unit ON (|).
- 4. Turn on power to the system control unit by toggling the power switch on the back of the unit ON (|).
- 5. Next, turn on the system control unit by pressing the power button on the front of the unit. This power button will glow when the unit is on.

The software will initialize and run automatically. The UI will indicate the software is INITIALIZING while the system checks for and establishes a connection with the attached peripheral cards. After the screen shows READY the system is ready for operation.

4.2.2 Controlling System Temperature

The PLATFORM temperature channel is used to drive control operations across the Cryostation, compressor, and vacuum control unit.

Cooling Down the System

- 1. In the UI for the CRYOSTATION instrument, navigate to the SAMPLE CHAMBER display screen and select the PLATFORM temperature channel to bring up its operation controls.
- 2. In the PLATFORM TARGET input box, enter the target temperature value. To reach the lowest possible base temperature, enter 0.0 K. Press SET to confirm.
- 3. Press the COOLDOWN command button. On the popup, confirm or adjust the custom parameters for the cooldown, then press COOLDOWN again to start.

The cooldown cycle will begin.

» NOTE

- The Stage 1 and Stage 2 temperatures will drop quicker than the platform temperature. The platform will drop faster after Stage 1 reaches 30 K.
- When the platform has reached a stable temperature, the ring surrounding the platform status readouts will change from flashing to solid.

Vacuum States During a Cooldown

- When a cooldown is initiated, the system will first establish a rough vacuum.
 - During this process, the system will automatically check for leaks. If a vacuum cannot be established, an error message will show in the UI and the cooldown process will be aborted. See System Diagnostics for troubleshooting information.
- After a proper vacuum level is established, the compressor will turn on; this may take up to 30 minutes.
- The pressure will continue to drop to a high vacuum state leveraging cryopumping.

» NOTE

- To achieve the best vacuum levels, set the cooldown target temperature to the system base temperature and allow the system to stabilize there before controlling to higher temperatures. This temperature set point is necessary to take full advantage of cryopumping.
- The VC1110 pressure gauge can only read out to ~0.1 mTorr. If pressure is below this value, the UI readout will indicate HIGH VAC.

Adjusting Platform Temperature

The platform temperature may be changed at any time by setting a new value using the PLATFORM TARGET input box. If the user is setting incremental temperature set points, the COOLDOWN and WARMUP buttons do not need to be pushed.

Warming Up the System

A warmup operation is used to bring the system to room temperature to access the sample chamber. The system may be warmed up actively with heaters at any time, including during a cooldown.

- 1. In the UI, navigate to the SAMPLE CHAMBER display screen and select the PLATFORM temperature channel to bring up its operation controls.
- 2. Press the WARMUP command button. On the popup, press WARMUP again to confirm.

An active warmup is much faster than using the STOP command. The heaters will automatically shut off when the entire system reaches 295 K, but the sample chamber will remain under vacuum.

Stopping a Cooldown or Warmup

At any time during a cooldown or warmup, the process can be stopped.

- 1. In the UI, navigate to the SAMPLE CHAMBER display screen and select the PLATFORM temperature channel to bring up its operation controls.
- 2. Press the STOP command button. On the popup, press STOP again to confirm.

This action will stop running the compressor/cryocooler and turn off any heater power going to the cryocooler stages and sample chamber platform. The system will begin to warm up naturally, but it will remain under a vacuum.

4.2.3 Venting and Pulling Vacuum

Vent to Access the Sample Chamber

After a cooldown and warmup cycle, the system remains under vacuum until the user is ready to access the sample chamber. To access the chamber:

- 1. In the UI, navigate to the SAMPLE CHAMBER display screen and select the PLATFORM temperature channel to bring up its operation controls.
- 2. Press the VENT command button. On the popup, confirm or adjust the custom parameters for the vent procedure, then press VENT again to confirm.

The sample chamber will be vented to atmospheric conditions.

» NOTE

Choosing to "Vent Continuously" will cause nitrogen (when attached) to flow through the chamber, even after the chamber is opened. This can help to keep the vacuum space as clean as possible during quick sample exchanges. To stop the flow of nitrogen when "Vent Continuously" is used, press the STOP command.

Keep the Sample Chamber Under Vacuum

When the system is not in use, it is recommended to keep the sample chamber under a vacuum to prevent moisture and contaminants from entering the sample space. To keep the system under vacuum:

- 1. In the UI, navigate to the SAMPLE CHAMBER display screen and select the PLATFORM temperature channel to bring up its operation controls.
- 2. Press the PULL VACUUM command button. On the popup, confirm or adjust the custom vacuum parameters, then press PULL VACUUM again to start.

The sample chamber will pull vacuum until the target vacuum pressure threshold is met.

» NOTE

Since this is an independent operation without cryopumping, the system will likely only be able to achieve a rough vacuum state. To stop the vacuum pump(s), press the STOP command. The sample chamber will remain at the current vacuum level until a COOLDOWN or VENT procedure is initiated.

4.3 Moving the System

If the system ever needs to be moved to a different lab or location, follow the steps below or reach out to the Montana Instruments Service team for guidance:

- 1. Save any important data on the system.
- 2. Power down the unit:
 - a. Tap (press and release) the power button on the front of the system control unit -or-
 - b. On the touchscreen UI, navigate to MENU > SYSTEM SETTINGS and select POWER OFF
- 3. Remove the helium and vacuum hoses, paying special attention to ensure the O-rings remain in place.
- 4. Remove the remaining cables and electrical connections.

If the system needs to be shipped, follow the remaining steps below

- 5. Reverse the steps in the <u>Installation Procedure</u>. Replace the red locking rings and c-spacers and lock the red shipping supports down on the tri-flange. Then unbolt the system from the optical table.
- 6. Re-pack the system in original packaging. Some components **must be** packed upright on a pallet to avoid damage.

NOTICE

- Follow all handling instructions for the individual components as outlined in the *General Hazard Information* section.
- If you are uncomfortable with moving the system on your own, or if you need to order any replacement packaging, please contact an authorized service representative.
- Do NOT attempt to disassemble any components of the system beyond the original state as shipped from the Montana Instruments factory.

Section 5 - Diagnostics & Troubleshooting

This section contains information for basic system diagnostics and troubleshooting advice. Diagnostics or repairs outside of the scope of this section should be completed by an authorized service representative.

5.1 Expected System Performance

Each system ships with a unique Certificate of Performance to demonstrate the factory performance of that Cryostation. Note that specifications and certifications represent factory acceptance testing performed on a bare platform without additional accessories or sample mounts.

The vibrations are measured in this configuration with a capacitive sensor that is also bolted to the table. The capacitive sensor measures the peak-to-peak vibrations in nanometers to a test fixture bolted to the sample platform.

5.2 System Diagnostics

5.2.1 Performance Issues

If a degradation in performance or other failures are experienced, check for these common issues:

Problem/Symptom	Possible Cause	Solution/Suggestion	
The system is unable to reach target temperature or cooldown takes longer than expected	See Temperature Optimization in section 5.3.1		
The system does not meet vibration specifications on the sample stage	See Vibration Mitigation in section 5.3.2		
The system will not pull rough vacuum – leak check failed	Vacuum leak	See Vacuum Check in section 5.3.3	
System condensing moisture on the sample, windows, exterior of			
the sample chamber, or vacuum case (black cylinder surrounding cryocooler). Exterior surfaces are cold to the touch.	Radiation (inner) and vacuum (outer) window are touching	Adjust windows so they do not touch. For low working distance setups, the windows may touch under vacuum, so be sure to check again after pulling the vacuum.	
Thermometer not working	Not properly installed or activated.	Make sure the temperature channel is enabled. Navigate to MENU >	
Heater not working	Not properly installed or activated.	INSTRUMENT SETTINGS > USER X. Toggle the Temperature Channel Enabled switch to the right to enable. Contact Customer Service for further troubleshooting & documentation	

5.2.2 Power and Communication Issues

If the system will not turn on, run commands, or display readouts, check for these common issues:

Problem/Symptom	Possible Cause	Solution/Suggestion
The system throws an electrical breaker	Wall power issue	1. Check the wall voltage using an AC voltmeter (digital multi-meter on the AC setting) to ensure it is in the defined ranges for each piece of equipment.
UI screen is black / does not turn on	Communication issue with system control unit	 Make sure all cables and power cords are connected properly. See Connecting System Cables and Power section of the Installation Procedure for details. Ensure the power switch on the back of the system control unit is ON (). Ensure the power button on the front panel is ON (glowing). Power cycle the system control unit. Toggle the power switch on the back of the unit OFF (o) then back ON (). Press the RESET button on the back of the UI touchscreen display.
	Touchscreen is faulty	 Contact MI Service for replacement Connect the external monitor to HDMI and connect a USB mouse to control the system OR Use VNC to control the system remotely
UI controls are frozen or non-responsive	Touchscreen display issue	1. Press the RESET button on the back of the UI touchscreen display.
The system does not initialize	Communication issue with the compressor	 Make sure all cables and power cords are connected properly. See Connecting System Cables and Power section of the Installation Procedure for details. Ensure the power switch on the back of the compressor is ON (). Ensure the ENABLE switch on the front panel is ON (). Power cycle the compressor. Toggle the power switch on the back of the unit OFF (o) then back ON ().

Vacuum pressure reading in UI indicates "SensorErr"	Communication issue with vacuum control unit		Make sure all cables and power cords are connected properly. See Connecting System Cables and Power section of the Installation Procedure for details. Ensure the power switch on the back of the vacuum control unit is ON (). Power cycle the vacuum control unit. Toggle the power switch on the back of the unit OFF (o) then back ON ().
---	--	--	--

5.2.3 Compressor Issues & Fault Messages

The variable-speed compressor has built-in mechanisms to check for issues. If the compressor alarm sounds or reads a fault message, check for these common issues:

For each of the errors below, the compressor may need to be restarted. To cycle the power:

- 1. Toggle the power switch on the back of the unit OFF (o) then back ON (|).
- 2. Ensure the ENABLE switch on the front panel is ON (|).
- 3. Ensure the helium pressure stabilizes in an acceptable range before proceeding (See *Helium Check section*).

Problem/Symptom	Possible Cause	Solution/Suggestion	
UI indicates "Cryocooler Disconnected"	Power dropped out with compressor or communication lost	Check cable and power connections. See <i>Power and Communication Issues section</i> .	
Compressor display screen blank		Check the EVENT LOG in the UI MENU for any fault messages.	
	Display screen failure	Contact Service	
System Halted Pressure Imbalance	Helium hose not completely connected	Check the hoses for leaks. See Helium Hose Fittings section.	
Return Pressure Too Low	Occurs when the return pressure is below its minimum value of 0.35MPa. The likely cause for this is the cryocooler stopped running or the return line is detached.	Check the connection on the return line and ensure the cryocooler is running Check the supply and return balanced pressures (1.78MPa)	
Supply Pressure Too High	Occurs when the supply pressure exceeds 2.75MPa. It is caused by similar conditions as the Return Pressure Too Low fault – the cryocooler stopped running or, in this case, the supply line is detached. In general, the Return Pressure Too Low fault will occur before this fault.	Check the connection on the supply line and ensure the cryocooler is running.	
Delta Pressure Too Low	Occurs when the difference between the supply and return line pressures falls below 0.3MPa. A problem occurred in the cryocooler causing the supply and return lines to approach a common value.	Check that the supply and return lines are attached correctly and not switched. If the compressor was running and stopped it will need time to rebalance the pressures.	
Capsule Over Temperature	Occurs when the capsule temperature exceeds 85°C. This could indicate improper oil flow or helium flow. It could also be caused by vents on the compressor being blocked.	Check that air flow through the compressor is not being inhibited by any external object.	
Helium Over Temperature	Occurs when the Cool Helium temperature exceeds 50°C. Similar to the capsule over-temperature fault, this is likely due to insufficient air flow through the cooling system.	Check that the compressor vents are clear of debris or impedances and that the fan is running.	
Exhaust Over Temperature	Occurs when the exhaust temperature of the cooling system exceeds 55°C. This is most likely due to insufficient airflow through the cooling system.	Check that the compressor vents are clear of debris or impedances and that the fan is running.	

Low Oil Flow	Occurs when the oil flow ratio falls below 0.4. This alarm indicates that the flow of oil through the oil cooling coil is insufficient. Since the oil flow depends on the differential pressure of the system, this alarm is commonly caused by low differential pressure. The differential pressure across the cryocooler depends on both the speed of the cryocooler motor and the compressor speed. A very high cryocooler speed could cause a drop in the differential pressure as could a very slow compressor speed. However, this scenario is highly unlikely.	Check that the supply and return lines are properly attached, and there is nothing that is causing a bypass of the cryocooler.
Oil Sensor Failure	Occurs when there is a problem with the oil sensor electronics. This is usually due to the oil sensor line being unplugged. This alarm could also be caused by a short in the sensor circuit. If this is suspected, contact Service	Check that the oil sensor wire harness connector is attached to the board.
Oil Return Failure	Occurs when the oil level falls below 30% in 30 seconds. This is indicative of a problem in the oil return line or in the oil level sensor itself.	Give the system a chance to settle and try again. If the problem continues, contact Service.
Adsorber Full	Indicates the adsorber needs to be replaced (occurs after ~30,000 hours of operation). Before this fault occurs, there will audible alarms & messages on the display to indicate a replacement may be needed soon.	Contact Service to schedule this replacement.
High Oil Level	If the oil return system is failing and the oil level continues to increase, this fault will occur when the oil level fills the pump.	Cycle the compressor power and restart the system. Contact Service if the alarm continues.

5.3 System Checks

There are several basic checks users can do to help diagnose general problems and help the system achieve optimal performance. Please handle these checks carefully and thoroughly, neglecting any one of these may have an impact on the base temperature or vibration performance.

5.3.1 Temperature Optimization

Cryostation platforms are optimized to control heat loads reaching the sample. To ensure the lowest possible base temperatures, follow the best practices below.

- 1. Use proper thermal lagging techniques for any wiring (including coax) entering the sample space. Ensure thermal clamps are properly tightened. See *Thermal Lagging Techniques in section* 3.5.2 for details.
- 2. Avoid "touches" between the various stages (such as a platform component touching a Stage 2 component, or a Stage 2 component touching a Stage 1 component), as these become conductive paths for heat transfer.
 - a. Ensure wiring or cabling does not contact the inside of the radiation shield.
 - b. Ensure the radiation shield or radiation windows do not contact any part of the outer vacuum housing or windows. For low working distance setups, confirm that windows do not touch after the vacuum is pulled.
- 3. Always use the appropriate wire material and gauge for the application. See <u>Sample Chamber</u> <u>Wiring in section 3.5</u> for details.
 - a. Avoid using copper wire unless required. The electrical conductivity of phosphor bronze is typically sufficient for most applications.
- 4. Ensure the screws holding the platform to the support base and the screws securing the radiation shield are in place and tightened (5 in-lb.). Failure to tighten the screws will reduce the ability of the system to pull heat from the sample platform and may increase vibrations.
- 5. Ensure a thin layer of N-grease or another thermal grease is used between metal-to-metal interfaces for proper thermal connection.
- 6. Use inner "cold" windows or blanks on the radiation shield whenever possible. Without them, the added heat load can significantly increase the base temperature of the platform.
- 7. Check the helium charge pressure. If there is a helium leak, cooling performance will be hindered significantly. See *Helium Check in section 5.3.4* for details.
- 8. Ensure the User temperature channel heater is NOT on.

» NOTE

During a cooldown, the system will steadily ramp down in temperature to 4.2 K, then can take longer to reach and stabilize at the base temperature. Options and energy inputs (laser power) may impact cooldown times and cause slightly higher base temperatures.

Identifying Heat Loads

The expected temperatures and temperature differences between Stage 1, Stage 2, Platform, and Sample under various operating conditions can be used to identify the source of an unwanted heat load. The temperatures below are examples to illustrate the troubleshooting process and should only be used as a general reference. It is recommended to compare with actual data from a previously successful cooldown on your system as a more accurate reference guide.

The base temperature for a standard CryoAdvance-50 under normal operating conditions is depicted below. The gradient between the platform and sample will change depending on the type of sample mount and other options.

	Stage 1	Stage 2	Platform	Sample
Temperature	27 – 28 K	2.4 – 2.7 K	2.7 – 3.0 K	2.8 – 3.2 K
Temperature Gradient			+0.2 – 0.4 K	+0.2 – 0.5K

Heat Load Between Stage 1 and Stage 2

In the example below, Stage 2 has a higher-than-normal temperature, as does the Platform and Sample. However, Stage 1 is colder than normal. The temperature gradient between Stage 2 and the Platform and the Platform and the Sample is normal. This indicates that the heat load is located between Stage 1 and Stage 2.

	Stage 1	Stage 2	Platform	Sample
Temperature	26 K	5.0 K	5.3 K	5.7 K
Temperature Gradient			+0.3 K	+0.4 K

Possible Cause: Crosslink rod touching the sidewall. Contact an authorized service representative for instructions to perform a touch test.

Heat Load Between Platform and Sample

In the example below, the Stage 2, Platform, and Sample temperatures are high. However, the most critical issue is the large temperature gradient between the Platform and Sample. The heat load is most likely located in between the Platform and Sample, and since the Sample temperature is high, the temperatures of the Platform and Stage 2 are also being pulled up.

	Stage 1	Stage 2	Platform	Sample
Temperature	28 K	3.0 K	3.3 K	5.0 K
Temperature Gradient			+0.3 K	+1.7 K

Possible Causes: The most likely culprits are those described in the *Temperature Optimization in section* 5.3.1, such as wires touching the radiation shield, wires touching the sample mount directly after the thermal clamp, use of improper wires, loose screws, missing N-grease, or missing radiation windows.

Heat Load Between Stage 2 and Platform

In rare cases, a higher-than-normal gradient may exist between Stage 2 and the Platform. This is typically caused by loose screws underneath the platform, which should not be accessed by users. If this heat load is present, contact an authorized service representative.

Heat Load on Sample Thermometer

If the sample is reading higher than normal, but all other temperatures are in a normal range, the most likely cause is an improperly mounted or lagged sample thermometer. Refer to *Mounting a Thermometer in section 3.3*.

5.3.2 Vibration Mitigation

The Cryostation and sample chamber have several vibration-damping design features to reduce the effects of cryocooler mechanical vibrations on the sample platform. To ensure the lowest possible mechanical vibrations, follow the best practices below.

- 1. Rigidly bolt the system to the optical table using all available mounting locations in the baseplate.
 - a. Bolts should be tightened sequentially, starting from one end of the Cryostation and progressing toward the other end
 - i. e.g. start at the sample chamber and move toward the back panel
- 2. Ensure all screws inside the sample space are tightened down.
- 3. Ensure that the red shipping rings and spacers beneath the shipping rings have been removed. The red shipping supports at the back of the Cryostation should be locked in the upper position off the tri-flange. See *Removing the Shipping Supports* section of the <u>Installation Procedure</u> for details. Having any of these components installed can introduce vibrations on the order of several microns.
- 4. Ensure the helium hoses are not touching the cabling or the optical table. They should also not be pressed against a wall or another surface. There should be a gentle 180° bend in the hoses to avoid any lateral tugging on the Cryostation as helium runs through them.
- 5. Ensure the vacuum hose is not in contact with anything moving, as this hose is somewhat springy.
- 6. Vibrational performance may vary depending on the operating speed of the compressor/cryocooler. Vibrations are minimized when using a variable speed compressor running at 22 Hz/50 Hz, which is automatically used during stable conditions at platform temperatures <20 K.
- 7. Ensure that optics are rigidly mounted to the table. Optics could be vibrating due to the Cryostation on the table and mounting them should be carefully considered to reduce this effect.
- 8. Check temperature performance. An increase in Stage2 temperature and a decrease in Stage1 temperature could indicate a crosslink rod touch that could inherently degrade vibrational performance.

5.3.3 Vacuum Check

The UI displays a pressure reading for the attached vacuum gauge sensor. If a system leak check fails, or if condensation or freezing is observed on or inside the chamber, a leak may be present. Please check the following:

- 1. Check to ensure the vacuum gauge properly reads atmospheric pressure (~500-770 Torr) when the system is at room temperature and vented. If it is not, the gauge may not be working properly.
- 2. Check that the vacuum housing and lid are in place and properly seated. Ensure no wires are pinched between the O-rings.

- 3. Check the O-rings on 1) the sample chamber vacuum housing 2) the vacuum housing lid and 3) the vacuum housing outer "warm" windows. Ensure the O-rings have a thin layer of L-grease and are completely free of debris or fibers.
 - a. All side panels also have an O-ring interface. Do NOT remove the side panels before consulting with an authorized service representative.
- 4. For the CryoAdvance-50 model, ensure the black Delrin screws holding the sample PCB in place are loose. If overtightened, these can cause the board to flex and compromise the O-ring seal.
- 5. Although nitrogen is optional, it helps keep the charcoal adsorbers clean and the inside of the Cryostation free from moisture. Using nitrogen is particularly important in humid environments.
 - a. Nitrogen purge cycles, as well as a platform bakeout, will help rid the chamber of contaminants before cooldown.
- 6. If a leak detector is accessible, use it to leak check the sample housing, Cryostation, vacuum hose, vacuum connections to find the source of the leak.

If the leak persists, please contact an authorized service representative.

There is enough charcoal in the system to freeze particles that may be introduced from a small leak. If there has been a large leak, or if a small leak persists over a period of time, the charcoal adsorbers will need to be recharged. To do this, run a COOLDOWN or PULL VACUUM operation with PLATFORM BAKEOUT (350 K for 60+ mins) and DRY NITROGEN PURGE (3+ times) enabled.

5.3.4 Helium Check

The helium pressure in the system can be checked when the compressor is idle.

- 1. Warm the system to room temperature and leave idle for at least one hour. Do not start a cooldown.
- 2. Ensure the power on the back of the unit is ON (|) and the ENABLE switch on the front panel is ON (|).
- 3. In the UI, navigate to the OVERVIEW display screen, then press the CRYOCOOLER reading to show compressor details.
 - a. Or, on the front panel of the compressor, toggle the DISPLAY button until it shows the supply and return values.
- 4. The values should be between 1.55 1.80 MPa (target 1.78 MPa) on both the supply and return. If the supply or return values are <1.55 MPa, helium will need to be recharged. See *Helium Recharge Process in section 5.3.4* for details.

Helium Hose Fittings


Improper connection or missing O-rings on the helium hose fittings can cause a loss in helium pressure and hinder the cooling performance.

To inspect fittings on the back of the compressor and Cryostation:

• Ensure the fittings are not cross-threaded.

- Ensure there is a **single** O-ring at each end of the hose and each connection point. The O-rings tend to dislodge from the hose and stay on the fitting (or vice versa).
 - o If this happens, the errant O-ring must be carefully removed and replaced in the correct location before reconnecting, otherwise it will not seat properly.



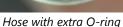



Extra O-ring removed



Left: Fitting missing O-ring Right: proper O-ring fitment

Helium Recharge Process

If the helium pressure is low (below 1.55 – 1.80 MPa), it should be recharged using a 99.999% (UHP) tank of pressurized helium. The accessory kit comes equipped with vent and recharge valves that can be used to flush the system and recharge. Contact an authorized service representative for instructions on this process, or find the service document "CryoAdvance B-01 Compressor Recharge Procedure" on the appropriate product page on the Montana Instruments website.

NOTICE

Follow instructions carefully

To avoid possible contamination, these instructions must be followed carefully.

A CAUTION

Ear protection recommended

Venting the helium lines typically creates a loud noise as pressure is released. Be prepared for venting noises when attaching the purge adapter to the helium lines.

Section 6 - Appendices

6.1 Related Documentation

For a copy of associated documentation, see below:

Document Title	Location
CryoAdvance-50 Documentation	https://www.montanainstruments.com/products/ca50r1
CryoAdvance-100 Documentation	https://www.montanainstruments.com/products/ca100r1
General Terms and Conditions of Sale	Direct Link to latest PDF Or find the link at the bottom of our homepage
Limited Warranty Agreement	https://www.montanainstruments.com/support/warranty-information
End-User License Agreement	Direct link to latest PDF