



BeatBox® Bead Remover - Quick Start Manual


This tool comprises the following items:

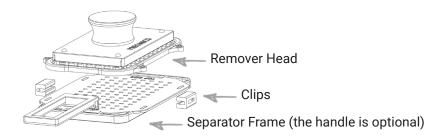

In addition, you will need the following items:

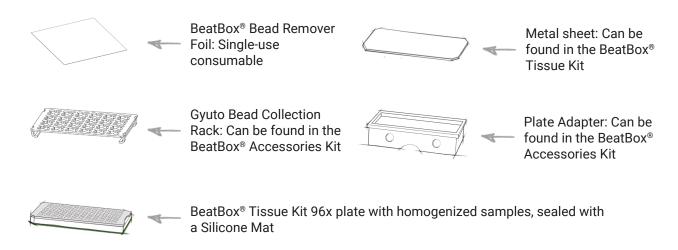

Prepare the plate

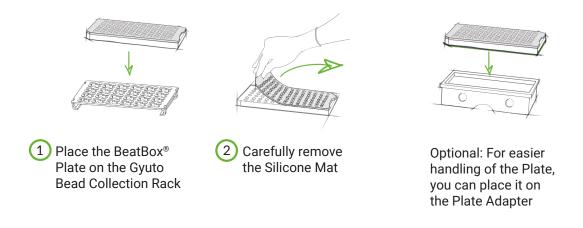

PREOMICS


Assemble the BeatBox Bead Remover




3 Optional: Attach the handle to the Separator Frame



Take the Separator
Frame, place a Foil on
the bottom side and
secure it with the clips.
Make sure that the part
within the outer frame is
covered by the Foil

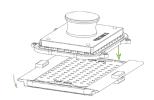


5 Turn the Frame around and place the Remover Head on top of it

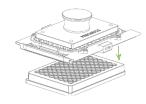

6 Place the BeatBox®
Bead Remover on top
of the BeatBox® Plate

Remove the Gyuto Beads

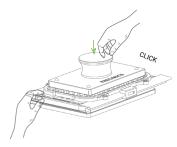


7 Hold the Separator Frame with one hand and press on the knob (standing position of the operator is recommended at this point)



8 Lift the tool off the plate by grabbing the knob with one hand and holding the plate in place with the other hand. The Gyuto Beads should sit on the Foil

Discard the Gyuto Beads

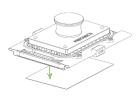

Place the BeatBox® Bead Remover with the Gyuto beads over the Metal Sheet



10 Separate the Remover Head from the Separator Frame by holding onto one of the clips and pulling on the round knob with the other hand



11 The Gyuto beads will stick to the Metal Sheet. Dispose of the Beads, Metal Sheet, and Foil according to regulations